

780 Turn On/Off Procedure

Kim Peifley

12/18/15

Procedure to Turn on LSM780

1. Turn on Main Switch.
2. Turn on System/PC Switch.
3. Turn on Computer.

1



2




3

Procedure to Turn on LSM780


4. Click LSM User to access desk top. No log in necessary. Wait until Windows fully opens before moving to step 5.
5. Turn on Components Switch.

4

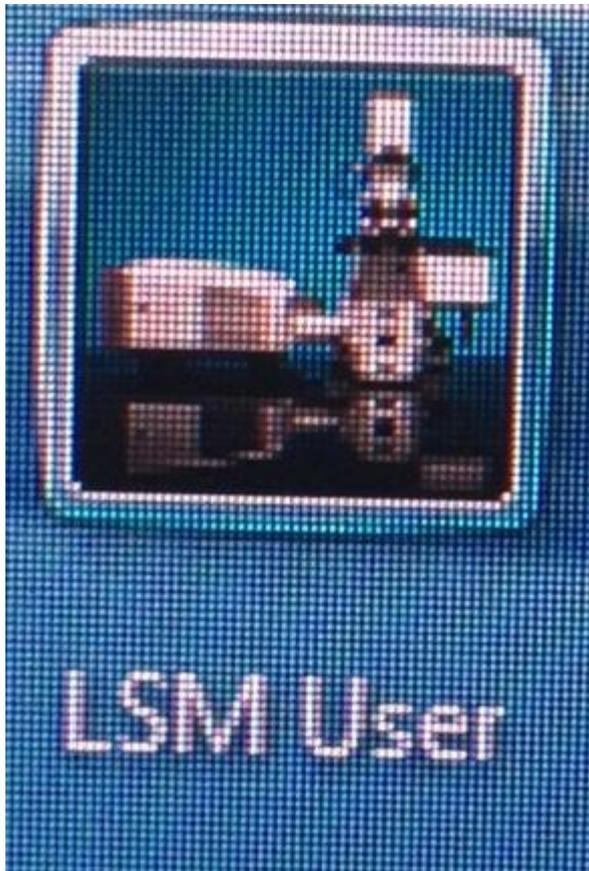

Procedure to Turn on LSM780

6. Turn on the X-Cite Lamp.
7. Turn Key to turn on Argon Laser Controller. Wait 5 minutes to warm up.
8. Flip toggle switch. Wait a few minutes until you see the light come on.
9. Adjust knob until it is at the point where red turns to green.

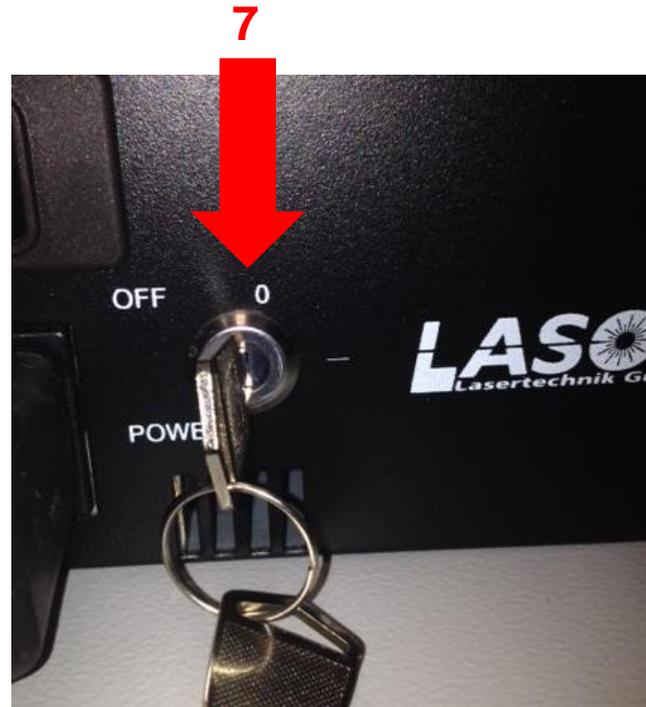

Procedure to Turn on LSM780


10. Click ZEN to open acquisition software.

10




Procedure to Turn Off LSM780

1. Close the software.
2. Move data to OMAL server.
3. Flip toggle switch for Argon Laser (see right).
4. Turn Key to turn off the Argon Laser Controller (see below)
5. Turn off Excite Lamp (see below).
6. Turn off Computer.

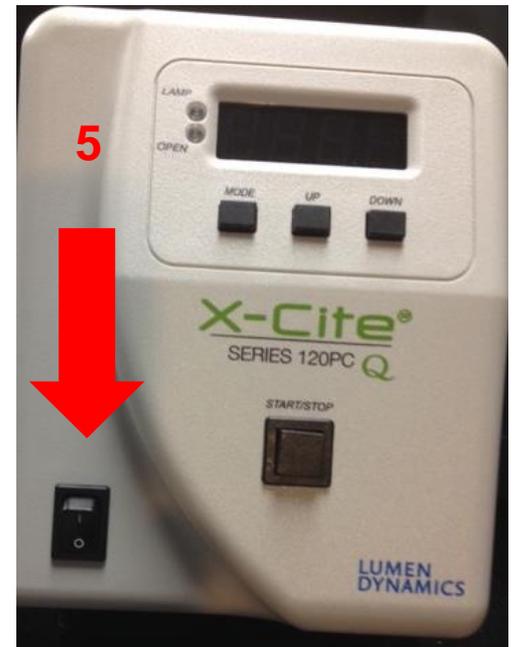

Procedure to Turn Off LSM780

7. Wait until you hear the fan shut off. This is cooling the lasers.
8. Turn off System/PC Switch (see below).
9. Turn off Components Switch (see below).
10. Turn off Main Switch (see below).

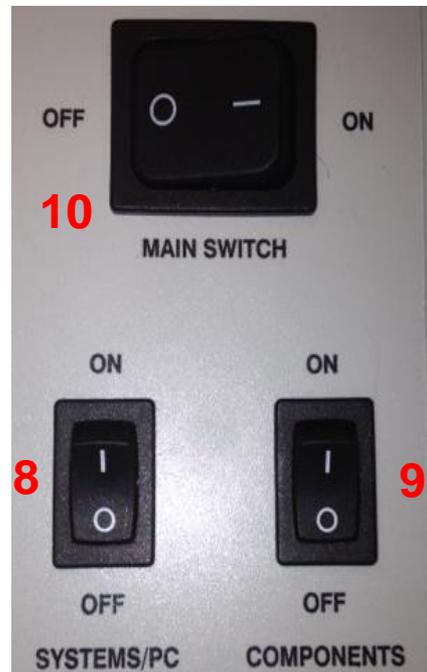

Troubleshooting

1. Instructions for when the Zen software on the 780 does not start up at all. In other words you only see a small window flash up for less than a second. This problem can happen if the system crashes when someone is using it (for example power outage). It is necessary to delete the folder illustrated in attachment.

